

MICROLIT PRO TIP™ | Micropipette Tips & Tip Boxes



ISO 17025



ISO 9001



ISO 13485

Experience Pipetting Precision with Microlit's ProTip™

Instructions for Use

Before using the tips, familiarize yourself with the operating manual of the relevant instrument.

Check the pack and contents carefully for any damage.

For filter tips, check whether there is a dual-filter in every tip and whether it is correctly positioned, undamaged and not twisted. If you detect any fault, do not use the tip.

The tips should always be emptied into a container so that no fluid sprays into the environment. After use, dispose of tips in accordance with relevant guidelines.

Tips should not be filled with fluids which attack Polypropylene (PP) or in the case of filter tips Polyethylene (PE) also.

If you are in any doubt as to the suitability of your product for its intended use or as to the reagent, you will find our contact details on the back page.



Manufacturer Recommended
Excellent compatibility with Microlit RBO & Microlit NERO - Assured accuracy and precision.



Exceptionally Clean Release
Considerably more hydrophobic than standard tips enabling even viscous fluids to be dispensed completely.



Certified free release
Free of detectable RNase, DNase, DNA, PCR inhibitors and tested non-pyrogenic.



Comprehensive Range
Both MICRO (up to 1000ul) and MACRO (up to 10ml) in standard (non filter, non sterile) and filtered varieties.



Flexible Packaging
Available in racked, reloads and bulk packaging.

SPECIFICATIONS & ORDERING INFORMATION

Product Description	Capacity	Pack/Box	Model No.
Micropipette Tip Packets (The tips are DNase, RNase and Pyrogen Free)	Upto 10 µl	100 tips/pack	MT/10/100
	Upto 200 µl	1000 tips/pack	MT/200/1000
	Upto 300 µl	1000 tips/pack	MT/300/1000
	Upto 1000 µl	1000 tips/pack	MT/1000/1000
	Upto 5000 µl	100 tips/pack	MT/5000/100
	Upto 10000 µl	100 tips/pack	MT/10000/100
Micropipette Tip Boxes (The tips are DNase, RNase and Pyrogen Free)	Upto 10 µl	96 tips/box	TPB-MT/10/96
	Upto 200 µl	96 tips/box	TPB-MT/200/96
	Upto 300 µl	96 tips/box	TPB-MT/300/96
	Upto 1000 µl	96 tips/box	TPB-MT/1000/96
	Upto 5000 µl	18 tips/box	TPB-MT/5000/18
	Upto 10000 µl	18 tips/box	TPB-MT/10000/18
Micropipette Filter Tip Packets (The filter tips are DNase, RNase and Pyrogen Free)	Upto 10 µl	100 tips/pack	MFT/10/100
	Upto 20 µl	100 tips/pack	MFT/20/100
	Upto 100 µl	100 tips/pack	MFT/100/1000
	Upto 200 µl	1000 tips/pack	MFT/200/1000
	Upto 300 µl	1000 tips/pack	MFT/300/1000
	Upto 1000 µl	1000 tips/pack	MFT/1000/1000
Micropipette Filter Tip Boxes (The filter tips are sterile, DNase, RNase and Pyrogen Free)	Upto 10 µl	96 tips/box	TPB-MFT/10/96
	Upto 20 µl	96 tips/box	TPB-MFT/20/96
	Upto 100 µl	96 tips/box	TPB-MFT/100/96
	Upto 200 µl	96 tips/box	TPB-MFT/200/96
	Upto 300 µl	96 tips/box	TPB-MFT/300/96
	Upto 1000 µl	100 tips/box	TPB-MFT/1000/100



OUR REACH



YOUR SATISFACTION IS OUR PRIORITY

Microlit products follow GMP norms and are manufactured under comprehensive QA (Quality Assurance) and QC (Quality Control). However, if you are dissatisfied with the operation of any of our products, get in touch with us or call your nearest Microlit dealer for free replacement.

CONTACT US

Microlit India

629 Pakramau, Kursi Road, Lucknow 226026, India
Phone: +91 9918625629, Email: info@microlit.com
www.microlit.com

Microlit USA

33 Wood Avenue South, Suite 600, Iselin, NJ 08830, USA
Phone: +1 732 321 0852, Email: info-usa@microlit.com
www.microlit.us

Microlit LATAM

Business Place Ibirapuera, Alameda dos Jurupis,1005,
Indianópolis, São Paulo, Brasil
Phone: +55 11 93269-5932, Email: info-latam@microlit.com